

ACTIVATED CARBON SAMPLE KIT FOR EARLY DRUG DEVELOPMENT

Benefits

- ◆ Activated carbon grades for adsorption of colorants, metals and process impurities
- ◆ Purify and synthesize more lead compounds in less time
- ◆ Maximize yield of target compounds while selectively removing unwanted impurities
- ◆ Improve efficiency by having convenient benchtop sample kit close at-hand
- ◆ Six high purity, half liter samples offered free-of-charge

Challenge and solution

In early drug development, process chemists in Chemical Research & Development departments are responsible for developing and optimizing a synthetic scheme and pilot procedure to scale-up promising lead compounds for clinical trials. Process chemists strive to maximize yield of the lead compound, while simultaneously removing unwanted impurities, including: 1) light, medium and dark colors; 2) metals including palladium, platinum, rhodium and ruthenium; and 3) process-related contaminants. Process chemists need to define a chemical process that is safe, reproducible and cost-effective.

It is almost impossible to predict which activated carbon grade will work best at removing unwanted impurities for any given sample. For this reason, we created a sample kit (Figure 1) for the pharmaceutical market containing six high purity activated carbon grades with unique selectivity for removing a range of the most common impurities (Figure 2).

The convenient design is ideal for placement on a laboratory benchtop. For each new sample, we recommend that process chemists try one chemically activated carbon and one steam activated carbon to determine which type works best, then optimize the purification method further by trying other grades in the kit.



Figure 1. Activated carbon sample kit for the pharmaceutical market

For the removal of	Product	Type	Activation	pH	Removal of color compounds			Activated carbon characteristics	
					Light colors (yellow)	Medium colors (light brown)	Dark colors (dark brown)	Purity	Filtration
Light/medium colors	NORIT® A SUPRA EUR/USP	PAC	Steam	Neutral	●	◐	◐	●	◐
	DARCO® G 60	PAC	Steam	5-8	◐	◐	◐	●	◐
	NORIT C EXTRA USP	PAC	Steam	Neutral	◐	●	◐	◐	◐
Medium/dark colors	NORIT CGP SUPER	PAC	Chemical	3.5-6	◐	◐	●	◐	◐
	NORIT CN1	PAC	Chemical	5.5-6	◐	◐	◐	◐	●
Light/dark colors	NORIT ROX 0.8	Extrudate	Steam	Neutral	◐	◐	◐	◐	N/A

Figure 2. Impurities removal table

● Best ◐ Better ◑ Good

Given that process chemists start with PAC but usually migrate to powdered activated carbon filter discs (Figure 3), we make these activated carbon grades available to the leading filter disc manufacturers, including: 3M, Pall, Millipore, Carlson, Gusmer, ErtelAlsop and Eaton Begerow. So whether you wish to use PAC for its ease of use and low cost, or prefer the cleaner filter disc product format to eliminate costly cleaning and validation steps, we have you covered.



Figure 3. Filter discs

How to order

The best way to receive a sample kit is to schedule a meeting with one of our Regional Sales Managers. We will deliver a presentation on activated carbon technology, its use in pharmaceutical applications and the kits will be handed out. Upon request, the presentation can also be tailored to meet your specific needs.

We offer two sample kits depending upon where you are located:

Sample item name	For region	Warehouse location	Key difference
Pharma Kit KZ	Europe, Middle East, Africa	Klazienaveen, Netherlands	Contains NORIT® SX PLUS LC
Pharma Kit MS	Asia Pacific, North America, South America	Marshall, TX, USA	Contains DARCO® G 60

Customers can also order the sample kit directly from our website cabotcorp.com/pharmasamplekit

Sample kit contents

1. Six half liter samples of activated carbon (5 powders and 1 extrudate)
2. USB drive containing pertinent documentation:
 - ◆ Product datasheets
 - ◆ Safety datasheets
 - ◆ Testing procedures
 - ◆ Product infosheet: Activated Carbon - Pharmaceutical Applications

Our sales, technical service and customer service teams are prepared to serve customers around the world. Contact us at cabotcorp.com/activatedcarboncontact



NORTH AMERICA
 Cabot Norit Americas, Inc.
 3200 University Avenue
 Marshall, Texas 75670
 USA
T +1 903 923 1000
F +1 903 923 1035

EUROPE, MIDDLE EAST & AFRICA
 Cabot Norit Nederland B.V.
 Astronaut 34
 3824 MJ Amersfoort
 THE NETHERLANDS
T +31 33 4648911
F +31 33 4617429

ASIA PACIFIC
 Cabot China Ltd.
 558 Shuangbai Road
 Shanghai 201108
 CHINA
T +86 21 5175 8800
F +86 21 6434 5532

SOUTH AMERICA
 Cabot Brasil Industria e Comercio Ltda.
 Rua do Paraiso 148 - 5 andar
 Sao Paulo 04103-000
 BRAZIL
T +55 11 2144 6400
F +55 11 3253 0051

The data and conclusions contained herein are based on work believed to be reliable, however, Cabot cannot and does not guarantee that similar results and/or conclusions will be obtained by others. This information is provided as a convenience and for informational purposes only. No guarantee or warranty as to this information, or any product to which it relates, is given or implied. This information may contain inaccuracies, errors or omissions and CABOT DISCLAIMS ALL WARRANTIES EXPRESS OR IMPLIED, INCLUDING MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE AS TO (i) SUCH INFORMATION, (ii) ANY PRODUCT OR (iii) INTELLECTUAL PROPERTY INFRINGEMENT. In no event is Cabot responsible for, and Cabot does not accept and hereby disclaims liability for, any damages whatsoever in connection with the use of or reliance on this information or any product to which it relates.

The NORIT and DARCO names are registered trademarks of Cabot Corporation or its subsidiary.

©2017 Cabot Corporation. All rights reserved worldwide.